The ERWEKA Export Manager

for direct PDF and XML export





Overview

Introducing the ERWEKA Export Manager: Simple and Fast Data Export! With the ERWEKA Export Manager, users can seamlessly export test protocols from devices SVM II, TAR II, EasyCheck, MultiCheck 6 and ZT 720/30 via their network, effortlessly converting them into PDF and XML formats. The Windows software, once installed on any PC within your lab network, automatically transforms your test report data into PDF and XML files. Say goodbye to time-consuming manual scanning of printed reports.

But that's not all. Our user-friendly interface allows you to define your desired destination folder, whether it is a local drive or a designated network location. Integrated into the labs own IT network, security measures and data compliance protocols valid for the network are also valid for those folders, ensuring valuable data always remains safeguarded. No compromises on data security or integrity.

Each exported file is specifically named, combining the device type, serial number, and the date and time of data reception. This means users will never lose track of valuable data, maintaining an effortless overview for easy accessibility and analysis.

The ERWEKA Export Manager is a light-weight solution, catering to both new and existing equipment. With its remarkable ability to receive data from multiple devices simultaneously, users will save time, effort, and money and streamline their laboratory workflows.

Supported devices

The Export Manager software supports the following ERWEKA devices:

- ZT 72x (from Firmware-Version 1.20) •
- ZT 73x (from Firmware-Version 2.02)
- EasyCheck (from Firmware-Version 1.38)
- MultiCheck 6 (from Firmware-Version 1.38)
- SVM II (from Firmware-Version 1.15)
- TAR II (from Firmware-Version 1.01)

Status

Shows the status of the network connection.

- Red: Values for "Port number" and "Target folder" are missing or incorrect.
- Yellow: The values entered are correct.
- Green: Data reception has been started. The software is ready to receive data from the device.

IP address

Shows the IP address of the PC on which the Export Manager is running. It is determined automatically.

Port number

Sets the number of the port where data is received from the device.

Supported data formats

The software supports ASCII and XML data formats for receiving data, and PDF and XML data formats for exporting.



Quick and easy PDF / XML export



Numerous supported **ERWEKA devices**



USB stick with software & copy protection dongle



Start listening

Switches the Export Manager to the "Waiting" status, in which data can be received. The button is active only after you have entered correct values for "Port number" and "Target folder".



3



Contact

Are you curious and want to find out more? Head over to our website and download our product brochures, watch videos of our equipment in action or find the ERWEKA dealer of your country.

+49 6103 92426-999



Fax:

support@erweka.com

E-Mail: sales@erweka.com Phone: +49 6103 92426-200



spareparts@erweka.com



www.erweka.com

www.facebook.com/erweka.gmbh

ERWEKA GmbH

Pittlerstr. 45 63225 Langen Germany E-Mail: sales@erweka.com Phone: +49 6103 92426-200 Fax: +49 6103 92426-999

Technical specifications of products described are stated without warranty and subject to change at any time without further notice. v.1.0.6.23

