
HORIBA


Explore the future

Laser scattering particle size distribution analyzer

LA-300

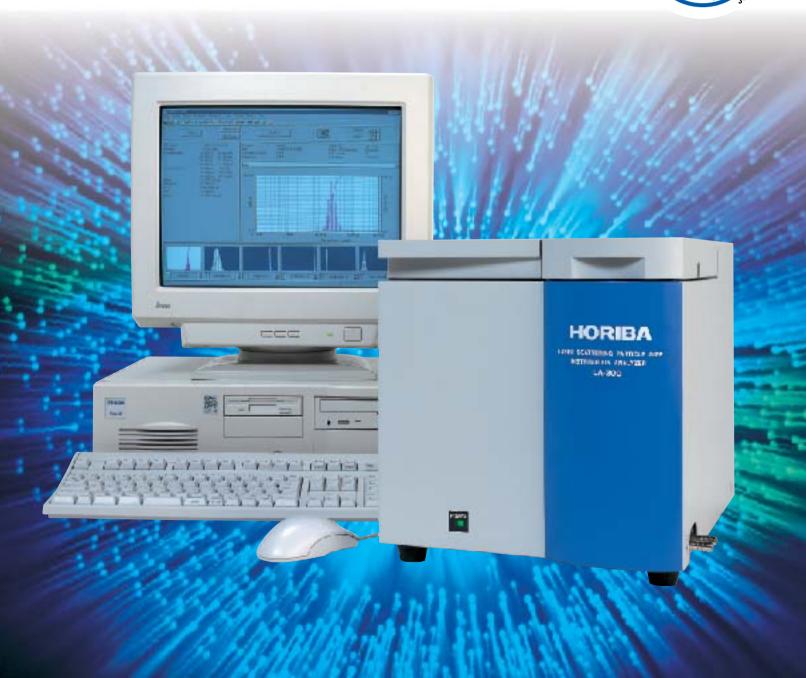



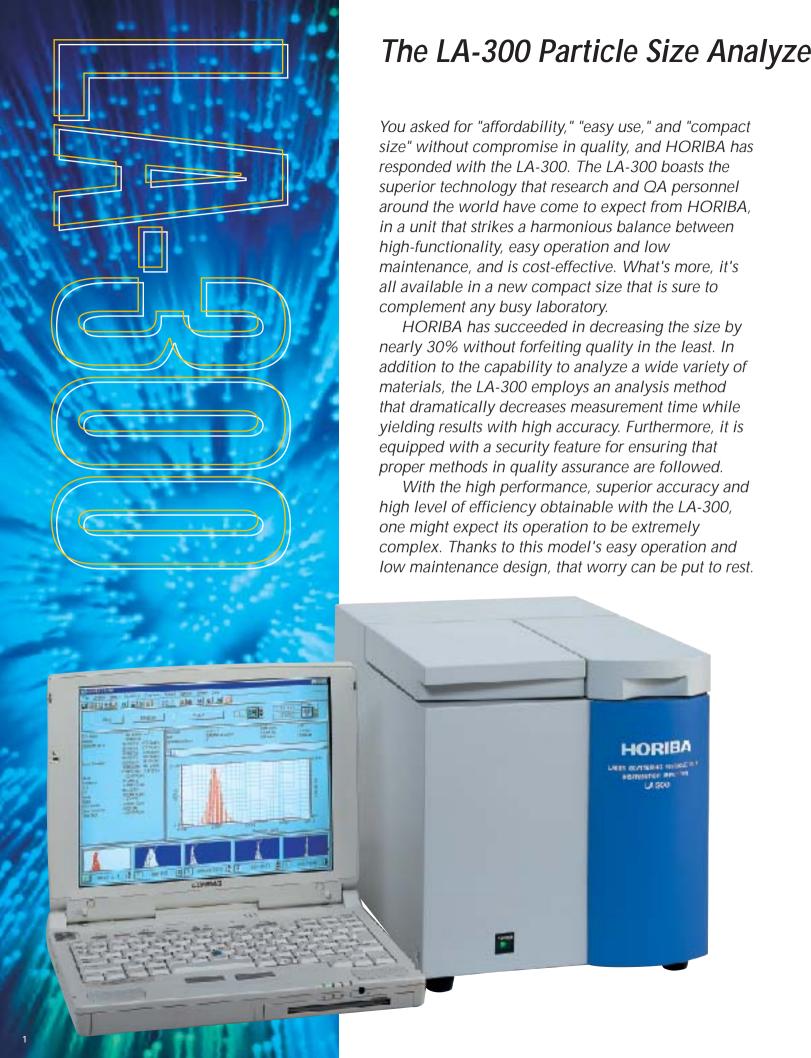


HORIBA's premium design in an economical package

High performance achieved in a compact size

The LA-300's revolutionary size makes it a welcome change to the other large and cumbersome designs, but rest assured that when it comes to performance, the LA-300 ranks right up there with the big guys.


A centrifugal pump system with an ultrasonic chamber makes it possible for measurements to be carried out at the same time as dispersion. Real-time displays of particle size distribution and sample concentration let the user check sample at all times. Maintenance is substantially reduced thanks to a strong centrifugal pump that resists residue buildup.

Using a diode laser, the LA-300 measures particles ranging anywhere between 0.1 and 600 μm in diameter with unprecedented precision, and processes high-resolution data for display within 20 seconds. The wide measurable range makes it ideal for application in fundamental studies and quality control of production lines in such diverse fields as pharmaceuticals, fine ceramics, metal powders, polymers, pigments, paints/coatings, adhesives, catalysts, minerals and clays, just to name a few.

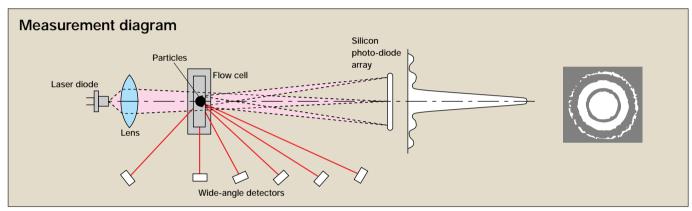

Measurements are performed in the LA-300 using a 650-nm laser diode (5mW). Six wide-angle detectors and a 36-channel ring-shaped silicon photo-diode array detector, which act as receptors for light refracting off of particles suspended in the flow cell, allow the LA-300 to process 42

data signals at once. Electrical signals corresponding to the intensity of the scattered light are used to calculate the size distribution of the particles. Based on the Mie scattering theory, this measurement method consistently yields superior repeatability with astounding precision.

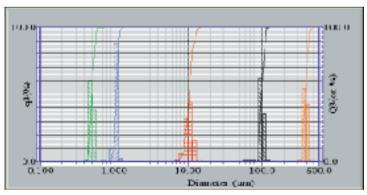

The above readouts were obtained from five separate analyses.


An example of the high-resolution readouts made possible with the 42-channel silicon photo-diode detector.

— HORIBA's premium design in an economical package

Designed for easy operation and maintenance

The software for the LA-300 offers extremely easy operation and low maintenance. A number of labor-saving features designed to facilitate operation are incorporated into the LA-300's application software.

A graphic user interface provides clear and concise dialog boxes, so performing operations requires little more than selecting the desired command buttons. Also included in this version are shortcut keys, convenient when performing main operations, and a window for quick assessment of memory availability.

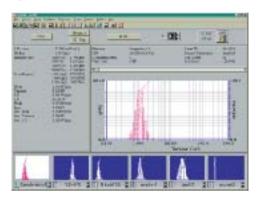

Auto-alignment

Thanks to a built-in auto-alignment function, the LA-300 is able to achieve high repeatability. This feature not only consistently maintains the ideal operation condition, but self-calibrates with the touch of a button, determining whether photo-diodes require re-positioning and adjusting them accordingly.

Multi-functionality at your fingertips

Security Feature


The LA-300's software application includes a window that allows the System Administrator to assign user-defined access for various LA-300 operations. The System Administrator can specify the attributes of user categories, as well as customize individual user IDs. User access is based on user name and password input, and can be limited to operations ranging from simply launching and closing the application, to more high-level functions, such as programming the LA-300's learning function, Learning Wizard.


Sequential Operation Wizard

Built into the LA-300 is HORIBA's unique operation function, Sequential Operation Wizard, which helps save time and labor when performing routine analyses. Once the user programs the desired measurement conditions and operation sequence into Sequential Operation Wizard, using an easy-to-follow graphic user interface, the recorded measurement sequence can be activated by just clicking on the appropriate button.

Print Layout

The Print Layout program in the LA-300 lets users perform custom layouts, complete with scaling and editing capabilities, for data output. Unlike most applications of this type, that offer a single printing format, the Print Layout application lets users edit the content, size and layout for displays of measurement conditions, graphs, results and other output data. Since a number of output formats can be saved, the format most appropriate for a particular print can be specified with ease. A preview screen is also provided so data can be viewed as it will be printed out.

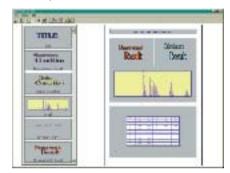

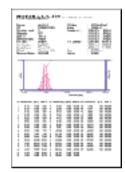

Low Maintenance

The LA-300's internal centrifugal circulation pump, for use with aqueous solutions, virtually eliminates the necessity for pump replacement since it is very difficult for powder residues to accumulate.

Additional features for meeting a variety of demands

Windows 98 Support

The LA-300 supports both Windows 95 and Windows 98. Now, data retrieved through a search program, for example, can be converted for processing in EXCEL and other Windows applications.

Flexible Data Processing

Various data processing options are provided to help you get the results you need. Among them is a graph display option that lets the user select to display a cumulative distribution graph, as well as selections for specifying histogram spacing, and how axes are to be displayed. Users can even choose to output to spreadsheets, or have graphs converted into bitmap or metafiles.

Production-boosting options



Autofill pump

This automatic fluid supply unit saves time and labor by automatically supplying the measuring unit with the sample fluid. It also provides automatic filling, rinsing of the fluid circulation system.



Fraction cell/cell holder


This unit is perfect for measuring very valuable or extremely small or hazardous samples. It allows liquid suspensions as small as 5 m ℓ to be measured.



For options that allow measurement of non-aqueous solutions, contact

HORIBA directly.

LA-300 with Autofill pump

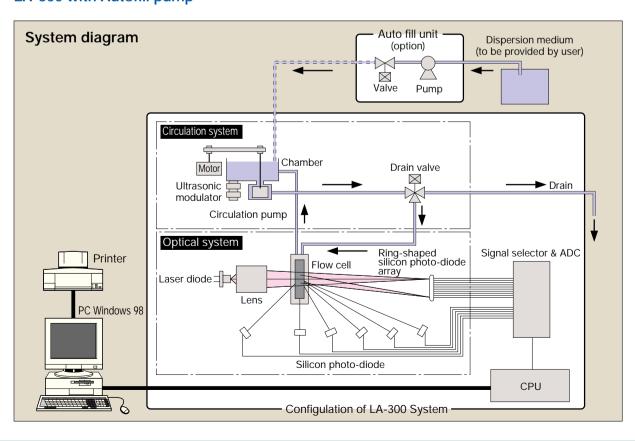

4

Specifications

LA-300 laser scattering particle size distribution analyzer

Principles of Measurement: Based on Mie scattering theory

Range of Particle Size Displayed: 0.1 to 600µm

Measurement Time:

Approximately 20 seconds, from start of measurement to display of data

Amount of sample required for measurement:

10 mg to 5 g

250 mL of dispersant aqueous* fluid when using the flow cell (differs with test samples)

Optical System:

Light source: 650-nm Laser diode, 5 mW

Photo-cell detector: 36-division ring shaped silicon photo-

diode forward detector array and 6 separate wide-angle silicon photo-

diodes

Sample Recirculation System:

Ultrasonic chamber, 15 W, 28 kHz

Centrifugal pump, discharge volume 5.5 L/min (in case of

distilled water), 15-step speed selection Sample flow cell, Tempax® glass

Power: 100/120/230 VAC ± 10%,

50/60 Hz, 150 VA

Communications: RS-232C

Proper Operation Temperature and Humidity:

15°C (59°F) to 35°C (95°F) less than 85% RH (no condensation)

External dimensions: 296 (W) x 420 (D) x 320 (H) mm

11.7 (W) x 16.5 (D) x 12.6 (H) in

Weight: Approximately 25 kg, 55 lb

Data processing, operating unit:

PC compatible computer running under PC/Windows 95 or

Windows 98

Printer

Monitor

Horiba continues contributing to the preservation of the global environment through analysis and measuring technology.



Options

Fraction cell and cell holder Cell material: Tempax® glass

Cell volume: 5 mL

Autofill pump


Flow rate: More than 1.2 L/min

Power: AC 100/120/230V ± 10%, 50/60 Hz, 80 VA

External dimensions:

165 (W) x 340 (D) x 320 (H) mm 6.5 (W) x 13.4 (D) x 12.6 (H) in Weight: Approximately 10 kg, 22 lb

External dimensions

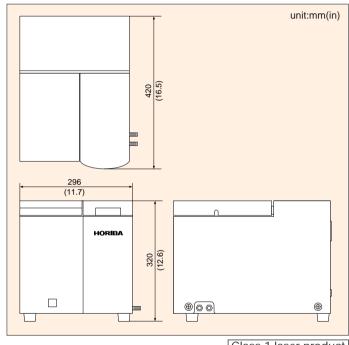

Class 1 laser product

* Contact HORIBA directly about options for use with non-aqueous solutions. Windows is a trademark of Microsoft Corporation



Please read the operation manual before using this product to assure safe and proper handling of the product.

- ●The contents of this catalog are subject to change without prior notice, and without any subsequent liability to this company.
- The color of the actual products may differ from the color pictured in this catalog due to printing limitations. • It is strictly forbidden to copy the content of this catalog in part or in full.

http://www.horiba.com e-mail: info@horiba.co.jp

●HORIBA, Ltd.— Head Office Miyanohigashi, Kisshoin

Minami-ku, Kyoto, Japan Phone: 81 (75) 313-8123 Fax: 81 (75) 321-5725

Tokyo Sales Office 1-7-8 Higashi-Kanda Chiyoda-ku, Tokyo, Japan Phone: 81 (3) 3861-8231 Fax: 81 (3) 3861-8259

Beijing Representative Office Suite 1409, Tower B, COFCO Plaza, No. 8, Jianguomennei Avenue. Beijing, China, 100005 Phone: 86 10-6522-7573 Fax: 86 10-6522-7582

Ann Arbor Facility 5900 Hines Drive Ann Arbor, MI 48108 U.S.A. INCORPORATED 1080 E. Duane, Suite. A Sunnyvale, CA 94086 U.S.A. Phone: 1 (734) 213-6555 Fax: 1 (734) 213-6525

Shanghai, China, 200020 Phone: 86 21-6415-3689/90 Fax: 86 21-6415-9746 ●HORIBA / STEC Kapianstrasse 5 A-3430 Tulin, Austria Phone: 43 (2272) 65225 Fax: 43 (2272) 65230

Phone: 1 (408) 730-4772 Fax: 1 (408) 730-8975

Shanghai Representative Office
Unit F1 16F
Unit F1 16F
No.15 Alley6, Lane 485,
Sec. 1, Kuang Fu Rd.,
No. 918, Huaihai Zhong Road,
Shanghai, China, 200020
Phone: 86 21-6415-3689/90
Fax: 886 (3) 5799164
Fax: 886 (3) 5799164

112-6 Sogong-Dong Choong-ku, Seoul, Korea Phone: 82 (2) 753-7911 Fax: 82 (2) 756-4972

●HORIBA KOREA Ltd.

● HORIBA INSTRUMENTS Pte. LTD.

10 Ubi Crescent #05-11/12, Ubi Techpark Singapore 408564 Phone: 65 6745-8300 Fax: 65 6745-8155

●HORIBA INSTRUMENTS LIMITED

Kyoto Close Summerhouse Road Moulton Park, Northampton NN3 6FL, U.K. Phone: 44 (1604) 542500 Fax: 44 (1604) 542699

●HORIBA INSTRUMENTS INCORPORATED

Irvine Facility 17671 Armstrong Avenue Irvine, CA 92614, U.S.A. Phone: 1 (949) 250-4811 Fax: 1 (949) 250-0924

●HORIBA EUROPE GmbH Head Office Hauptstrasse 108

D-65843 Sulzbach Germany Phone: 49 (6196) 6718-0 Fax: 49 (6196) 641198 Leichlingen Facility Julius-kronenberg Strasse D-42799 Leichlingen Germany Phone: 49 (2175) 8978-0 Fax: 49 (2175) 8978-50

HORIBA FRANCE Rue L. et A. Lumière Technoparc F-01630 St-Genis-Pouilly

France Phone: 33 (4) 50-42-27-63 Fax: 33 (4) 50-42-07-74

HORIBA SWEDEN

●HORIBA GmbH

Kaplanstrasse 5

Hertig Carlsvag 55-57 S-15138 Södertälje Sweden Phone: 46 (8) 550-80701 Fax: 46 (8) 550-80567

HORIBA CZECHIA Organizachi slozka Praha Petrohradska 13 CZ-101 00 Praha 10, Czech Republic Phone: 420 (2) 717-464-80 Fax: 420 (2) 717-470-64

HORIBA ITALY Europalace Corso Torino 43/45 10043 Orbassano, Torino, Italy Phone: 39 (011) 9040601 Fax: 39 (011) 9000448

Bulletin:HRE-3643C