

LM10 Microfluidizer™ High Shear Fluid Processor

Digitally Controlled Lab Unit For Small Sample Material Processing

Microfluidizer™ technology converts fluid pressure more efficiently into shear forces, leading industry performance standards in high pressure homogenization.

A unique solution to maintaining consistent process pressure, ensures 100% of your material gets exactly the same treatment. So whether you are working with small lab samples or production volumes, the Microfluidizer™ is unmatched in submicron size reduction/disruption, product yield, and guaranteed process scale-up.

Recommended For:


- Emulsions
- Dispersions
- Liposomes
- Cell Disruption



Model shown is subject to change depending on options selected

Unique Benefits of the LM10

- Achieve unmatched size reduction/disruption performance at lower process pressure:
 - Limit sample temperature rise during processing
- Higher product yields after sterile filtration with tighter particle size distribution
- Save on development time in pilot/production with linear volumetric scale-up
- Enhanced repeatability with easy-to-use digital pressure control
- Improved experimental reliability with real-time temperature monitoring
- Ensure dependable process performance over time with maintenance reminder and operator alerts
- Adapt to future experimentation possibilities with field upgradable firmware

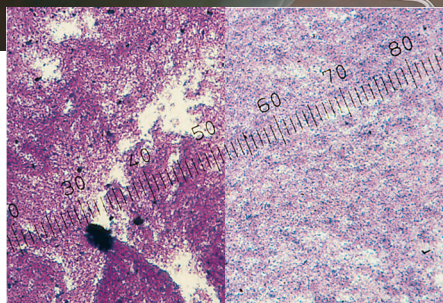

Before Processing After Processing

Along with other uses, the patented Microfluidizer™ can also be used for fine particle deagglomeration.



Operating Principle

The LM10 contains an intensifier pump designed to supply the desired pressure at a constant rate to the product stream. As the pump travels through its pressure stroke, it drives the product at constant pressure through precisely defined fixed-geometry microchannels within the interaction chamber.

As a result, the product stream accelerates to high velocities, creating shear rates within the product stream that are orders of magnitude greater than any other conventional means. All of the product experiences identical processing conditions, producing the desired results, including: uniform particle and droplet size reduction (often submicron), deagglomeration and high-yield cell disruption.

A removable cooling coil and a cooling bath are incorporated into the design to promote optimal temperature control.

Standard Features



- 300 ml glass reservoir
- Ceramic interaction chamber assembly
- Cooling coil and chilled bath assembly
- Thermocouple at product outlet
- Easily autoclavable

Options

- Diamond interaction chambers
- Process pressure transducer
- Larger capacity glass or stainless steel reservoirs
- Thermocouple at product inlet
- Process fluid recirculation assembly/kit
- Gasket materials

Specifications

Pressure Range	Up to 23,000 psi (1,546 bar)
Minimum Sample Size	14 ml
Flowrate Range	250 - 600 ml/min
Feed Temperature	Maximum of 165°F (75°C)
Electrical Power Requirement	85-260VAC, 50/60Hz, 5A
Pump Actuator	Pneumatically Driven
Air Compressor Requirement	57 scfm @ 120 psi (27 l/s @ 8.3 bar)
Noise Level	<80dB TWA
Dimensions	19"H x 19"W x 27"L (48cm x 48cm x 69cm)
Weight	75 lbs (34 kg)

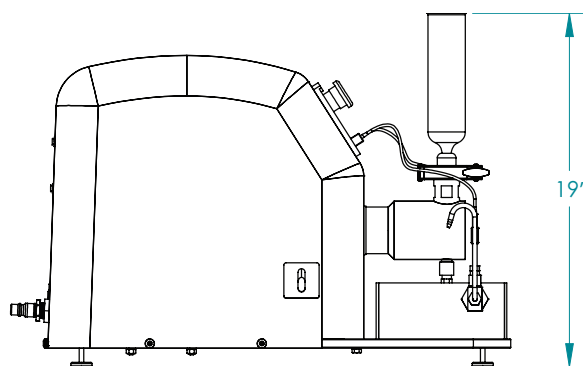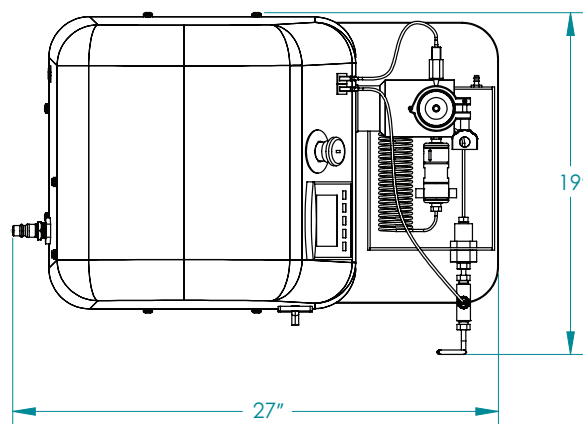

*Dimensions Are Approximate

Microfluidics

Microfluidics International Corporation
90 Glacier Drive, Suite 1000 • Westwood, MA 02090, USA
Tel: 617-969-5452 • 800-370-5452 • Fax: 617-965-1213
Email: mixinginfo@idexcorp.com • www.microfluidicscorp.com

Distributed in Canada by:
ATS Scientific Inc.
4030 Mainway
Burlington ON L7M 4B9
1-800-661-6700
www.ats-scientific.com



©2014 by Microfluidics. All rights reserved. 01/14 IH .5K